VEVOR[®]

TOUGH TOOLS, HALF PRICE

Technical Support and E-Warranty Certificate <https://www.vevor.com/support>

ELECTRONIC BALANCE

USER MANUAL

MODEL: HZ-B50002,HZ-B5002,HZ-B3002

We continue to be committed to provide you tools with competitive price.

"Save Half", "Half Price" or any other similar expressions used by us only represents an estimate of savings you might benefit from buying certain tools with us compared to the major top brands and does not necessarily mean to cover all categories of tools offered by us. You are kindly reminded to verify carefully when you are placing an order with us if you are actually saving half in comparison with the top major brands.

VEVOR®

TOUGH TOOLS, HALF PRICE

ELECTRONIC BALANCE

MODEL: HZ-B50002,HZ-B5002,HZ-B3002

NEED HELP? CONTACT US!

Have product questions? Need technical support? Please feel free to contact us:

 CustomerService@vevor.com

This is the original instruction, please read all manual instructions carefully before operating. VEVOR reserves a clear interpretation of our user manual. The appearance of the product shall be subject to the product you received. Please forgive us that we won't inform you again if there are any technology or software updates on our product.

 **WARNING:**

Please read this manual carefully before using the product. Failure to do so may result in serious injury. SAVE THIS MANUAL

Do not put in flammable liquid.

- Could cause a fire.

Do not spray water or liquid over the device which will causes electric shock . fire and malfunction.

Do not use the machine at any damp places.

Could cause electric shock fire and malfunction.

Do not touch the electrical connection with any metal objects.

- Could cause electric shock . fire and malfunction.

Do not disassemble , repair , modify the machine.

Could cause electric shock or injury.

Do not use products other than special adapter.

- could cause fire electric shock.

Do not touch the power plug with wet hands , especially when inserting or removing the plug.- Could cause fire . electric shock and malfunction.

Remove the power plug from the socket if the device is not going to be used for a long time.- Easy to cause fire , electric shock and malfunction.

Precautions for Use

OPERATION

1. Place the balance on a smooth and steady table. Please adjust balance through adjust screws at the bottom of the balance if the table is not smooth or steady. Avoid vibration, sunshine, air movement and magnet.
2. The air temperature should be kept at about 5C-35C at all times with a relative humidity level of about 50-85%.
3. The balance accepts AC input voltages of 220V, DC input voltages of 9V.
4. Please connect power supply and turn the balance on. It will display "8.8.8.8.8.8.", "Max. Capacity" and "_" step by step. You can use it after it displays "0.0", "0.00" or "0.000". (You had better warm the balance up for half an hour before using.)
5. Place the object you want to weigh on the balance pan, it will display the weight value.
6. Place a container on the balance pan if you want to weigh the object located in a container. Press "TARE/ZERO" key to zero the display. Place the object in the container, the balance will display the weight value.
7. Press "PCS" key for part count, the balance will display "PCS" and "10". Choose average number from "10", "20", "50", "100", "150", "200", "250" to "500" by pressing "CAL", Place "10", "20", "50", "100%", "150", "200", "250" or "500" pieces of same object you want to weigh on the balance pan, press "PCS" key to confirm. Please take the objects away from the balance pan. Now you can count object. Press "PCS" key if you want to escape count mode.
8. Press "UNIT" key for choosing weighing unit: g, ct, oz, lb.

LINEARITY CALIBRATION

Press and hold "CAL" for 5 seconds, the balance displays "CAL" and a weight value, Place relevant weights on the balance pan according to the weight values. At last, remove the weight, it displays "0.0", "0.00" or "0.000", You had finished the calibration.

SAVE THESE INSTRUCTIONS

TECHNICAL SPECIFICATIONS

Model	HZ-B50002
Type	Square scale pan,LCD display screen
Range(g)	5000 g
Accuracy(g)	0.01 g
Size(mm)	260*200*70
Net Weight (kg)	1.9
Gross Weight (kg)	2.2

Model	HZ-B5002
Type	Circular scale pan,LCD display screen
Range(g)	500 g
Accuracy(g)	0.01 g
Size(mm)	260*200*70
Net Weight (kg)	1.6
Gross Weight (kg)	2.0

Model	HZ-B30002
Type	Square scale pan,LCD display screen
Range(g)	3000 g
Accuracy(g)	0.01 g
Size(mm)	260*200*70
Net Weight (kg)	1.6
Gross Weight (kg)	2.3

*Products such as specifications , the appearance and the design are subject to modification without prior notice.

NOTICE:

1. The balance will display “_ - - ” if your object overloads 0.2 percent of the full capacity, please remove your object that time.
2. Please switch the balance pan according to the direction of hour hand if you want to remove the pan, Do not lift it up2directly.
3. Error Information Table.

Error Information	Specification	Transaction
_____	It means that the object's weight on the balance pan is less than 0 gram	Calibrate the balance
- - - - -	It means that the object's weight on the balance pan is overload	Calibrate the balance.
Err-1	It means that you turn on and off the balance continuously	Turn the balance off then turn it on after 3 seconds
Err-2	It means that the object's weight is not stable	Do not operate anything

4. For more details of technology, contact manufacturer please.

Address: Baoshanqu Shuangchenglu 803long 11hao 1602A-1609shi
Shanghai

Imported to AUS: SIHAO PTY LTD, 1 ROKEVA STREETEASTWOOD
NSW 2122 Australia

Imported to USA: Sanven Technology Ltd., Suite 250, 9166 Anaheim
Place, Rancho Cucamonga, CA 91730

Pooledas Group Ltd
Unit 5 Albert Edward House, The Pavilions
Preston, United Kingdom

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VEVOR[®]

TOUGH TOOLS, HALF PRICE

Support and E-Warranty Certificate

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Importado a AUS: SIHAO PTY LTD, 1 ROKEVA STREETEASTWOOD NSW 2122
EE. UU.: Sanven Technology Ltd., Suite 250, 9166 Anaheim Place, Rancho Cucamonga, CA 91730
Grupo Pooledas Ltd.
SHUNSHUN GmbH