VEVOR[®]

TOUGH TOOLS, HALF PRICE

Technical Support and E-Warranty Certificate

www.vevor.com/support

OVERHEAD MIXERS

USER MANUAL

MODEL NO.: 98-I-AB-1000, 98-I-AB-500, 98-II-AB-2000

We continue to be committed to provide you tools with competitive price.

"Save Half", "Half Price" or any other similar expressions used by us only represents an estimate of savings you might benefit from buying certain tools with us compared to the major top brands and does not necessarily mean to cover all categories of tools offered by us. You are kindly reminded to verify carefully when you are placing an order with us if you are actually saving half in comparison with the top major brands.

VEVOR®

TOUGH TOOLS, HALF PRICE

Magnetic Stirrers

MODEL NO.: 98-I-AB-1000, 98-I-AB-500, 98-II-AB-2000

NEED HELP? CONTACT US!

Have product questions? Need technical support? Please feel free to contact us:

Technical Support and E-Warranty Certificate
www.vevor.com/support

This is the original instruction, please read all manual instructions carefully before operating. VEVOR reserves a clear interpretation of our user manual. The appearance of the product shall be subject to the product you received. Please forgive us that we won't inform you again if there are any technology or software updates on our product.

WARNING:

Please read this manual carefully before using the product. Failure to do so may result in serious injury.

	Warning! ● Read the operating instructions carefully before use. ● Ensure that only trained staff works with the instrument.
	Protective ground contact! ● Make sure that socket is grounded (protective ground contact) before use.
	Risk of burn! ● Caution when touching the housing parts and the hotplate which can reach a temperature of exceed 300 °C. ● Pay attention to the residual heat after switching off.

- When operating wear personal safety protection to avoid the risk of:
 - Splashing and evaporation of liquids
 - Release of toxic or combustible gases
- Set up the instrument in a spacious area on a stable, clean, non-slip, dry and fireproof surface. Do not operate the instrument in explosive atmospheres, with hazardous substances or around water.
- Gradually increase the speed, reduce the speed if:
 - Stirring bar breaks away due to high speed
 - The instrument is not running smoothly, or the container is not centered on the plate
- Temperature must always be set to at least 50°C lower than the flash point of the media used.
- Be aware of hazards due to:
 - Flammable materials or media with a low boiling temperature
 - Overfilling of media
 - Unsafe container
- Process pathogenic materials only in closed vessels.
 - If the case of the stirrer bar is PTFE, please note:
 - Elemental fluorine, three fluoride and alkali metals will corrode the

PTFE and Halogen alkanes making it expand at room temperature.

- Molten alkali, alkaline earth metals or their solution, as well as the powder in the second and third ethnic of the Periodic Table of Elements will have a chemical reaction with PTFE when temperature reaches 300 ~ 400°C.
- Check the instrument and accessories prior to each use. Do not use damaged components. Safe operation is only guaranteed with the accessories listed in the “Accessories” section. Accessories must be securely attached to the device and must be removed when not in use. Always disconnect the power before fitting accessories.
- When the external temperature sensor is needed, the tip of the measuring sensor must be at least 5-10 mm from vessel bottom and wall.
- The instrument can only be fully disconnected from the main power supply by turning off the main or disconnecting the plug.
- The voltage stated on the label must correspond to the main power supply.
- Ensure that the main power supply cable does not touch the hotplate. Do not cover the device.
- The instrument must only be operated by experts.

Keep away from high magnetic fields.

SAVE THESE INSTRUCTION

PROPER USE

The instrument is designed for mixing and/or heating liquids in schools, laboratories or factories

- Observe the minimum distances between multiple units, and distances to the rear wall and above the assembly (min. 100 mm).

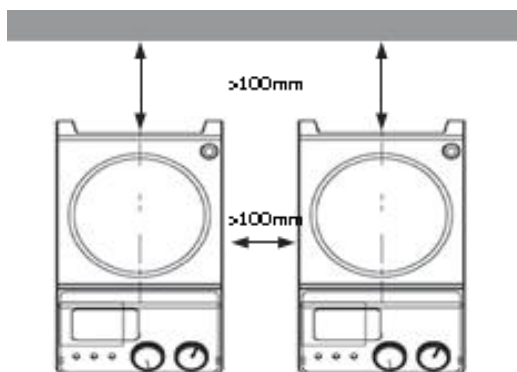
Figure 1

This device is not suitable for use in residential areas or other constraints mentioned in Safety Instructions.

TECHNICAL SPECIFICATIONS

Model	98-I-AB-1000	98-I-AB-500	98-II-AB-2000
Voltage [VAC]	220~240	220~240	220~240
Frequency [Hz]	50	50	50
Power [W]	350	250	500
Maximum stirring capacity [mL]	1000	500	2000
Maximum heating temperature [°C]	380	380	380
Product Dimensions (L*W*H)[mm]	φ285*210	φ220*165	φ330*230
Packing Size (L*W*H) [mm]	300*300*270	230*230*230	330*330*280
Net Weight [kg]	3.5	2	4.5
Gross Weight [kg]	4.2	2.5	5.5

*Products such as specifications, appearance, and design are subject to modification without prior notice.

PACKAGE CONTENTS

MODEL NO.: 98-I-AB-1000, 98-I-AB-500

Machine*1

Power cord*1

Stirrer bar*1

MODEL NO.: 98-II-AB-2000

Machine*1

Temperature
probe*1


Power cord*1
Stirrer bar*1

1 set of
brackets tools

PRODUCT FUNCTIONS


MODEL NO.: 98-I-AB-1000, 98-I-AB-500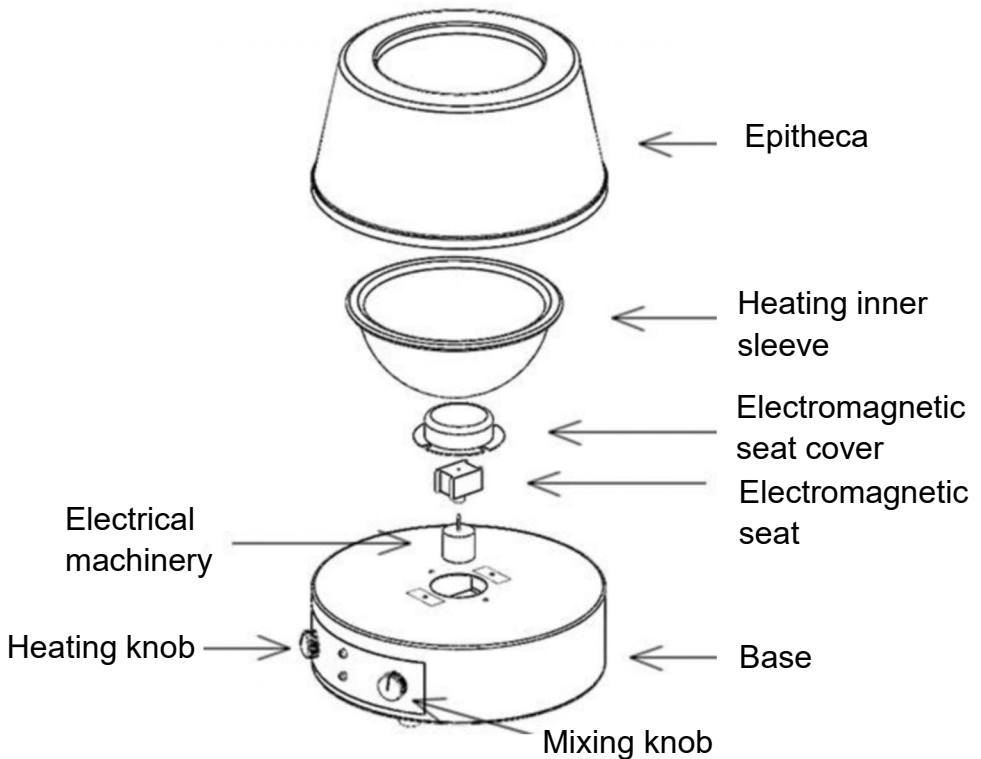

Operation mode:

1. Place the machine on a horizontal desktop, and then connect the power cord to the power outlet on the back of the base.
2. Plug the other end of the power cord into a standard socket.
3. Place the machine in a stable position, put the flask containing liquid medicine in the center of the working plate, put the appropriate rotor in the flask, and make it in the center of the flask,
4. turn on the power, adjust the temperature by rotating the "HEAT" knob at different angles, the "HEAT" indicator light is on when heating, and the heating power is adjusted by the "HEAT" knob.
5. Adjust the stirring speed by rotating the "STIR" knob at different angles, the "STIR" indicator light is on when stirring, and the STIR speed is adjusted by the "STIR" knob.


MODEL NO.: 98-II-AB-2000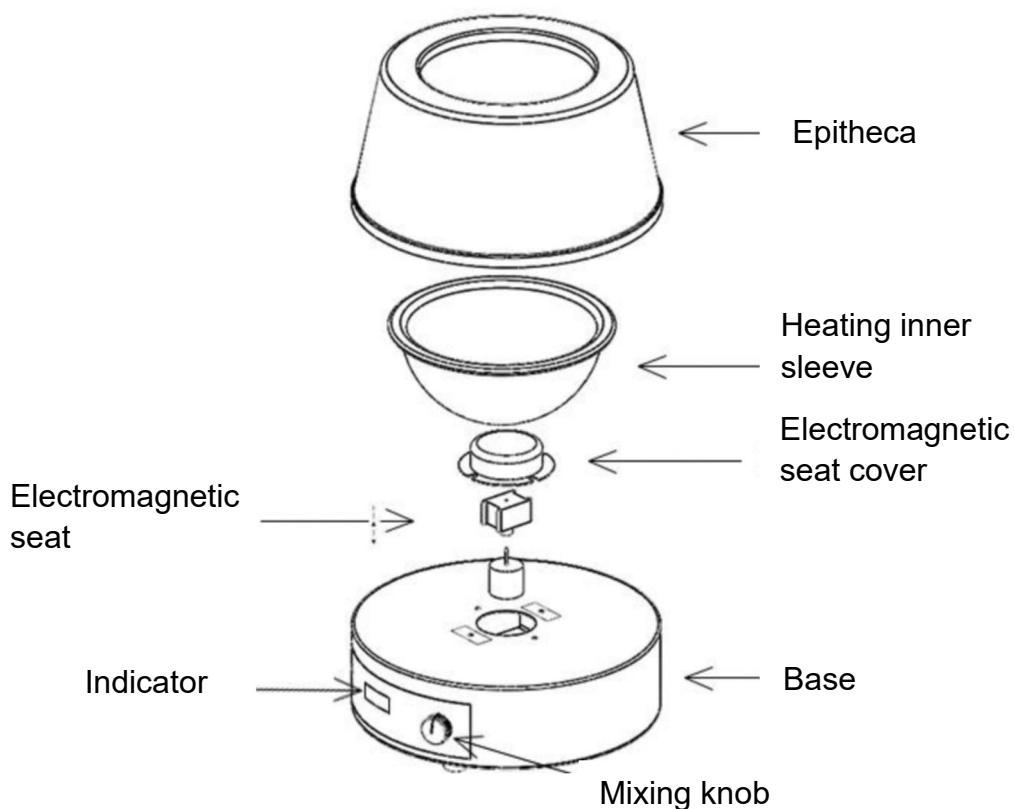

Operation mode:

1. Place the machine on a horizontal desktop, and then connect the power cord to the power outlet on the back of the base.
2. Plug the other end of the power cord into a standard socket.
3. Pay attention to the power switch on the rear side of the machine base. (-) means open, (o) means close. (There is no switch on Double-knob type Heating Mantle).

4. Place the machine in a stable position, put the flask containing liquid medicine in the center of the working plate, put the appropriate rotor in the flask, and make it in the center of the flask, turn on the power, press "S"key to enter the setting state, press "Increase" or " Decrease " key to adjust the temperature, and then the machine starts to heat. Turn the "STIR" knob to make the machine start to stir , and rotate the"STIR" knob at different angels to adjust the rotational speed.

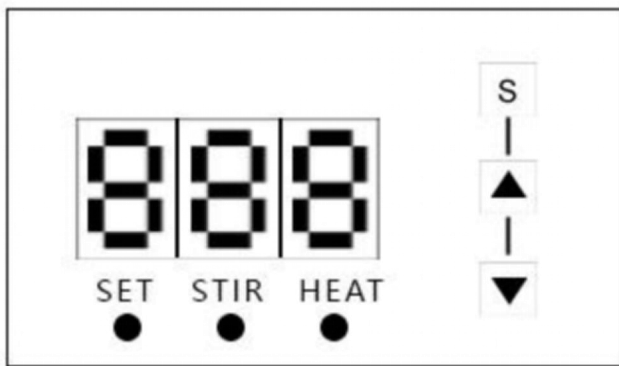5. If need to measure the temperature of the liquid in the flask, connect the bracket, insert the sensor lead plug into the sensor jack on the back of the machine, and then insert the sensor into the liquid, the temperature displayed on the meter at this time is the temperature of the liquid in the flask. (There is no external sensor for Double-knob type Heating Mantle).

This product is subject to the provision of European Directive 2012/19/EC. The symbol showing a wheeled bin crossed through indicates that the product requires separate refuse collection in the European Union. This applies to the product and all accessories marked with this symbol. Products marked as such may not be discarded with normal domestic waste, but must be taken to a collection point for recycling electrical and electronic devices.

Manufacturer: Shanghaimuxinmuyeyouxiangongsi
Address: Shuangchenglu 803nong11hao1602A-1609shi, baoshanqu, shanghai 200000 CN.
Imported to AUS: SIHAO PTY LTD. 1 ROKEVA STREETEASTWOOD NSW 2122 Australia
Imported to USA: Sanven Technology Ltd. Suite 250, 9166 Anaheim Place, Rancho Cucamonga, CA 91730

UK	REP
----	-----

YH CONSULTING LIMITED.
C/O YH Consulting Limited Office 147, Centurion House, London Road, Staines-upon-Thames, Surrey, TW18 4AX

EC	REP
----	-----

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MODELLO NUMERO: 98-II-AB-2000
Importato in AUS: SIHAO PTY LTD. 1 ROKEVA STREETEASTWOOD NSW 2122 Australia
Importato negli USA: Sanven Technology Ltd. Suite 250, 9166 Anaheim Place, Rancho Cucamonga, CA 91730
E-CrossStu GmbH