VEVOR

Affordable. Reliable. Home Improvement.

FLAMMABLE CABINET

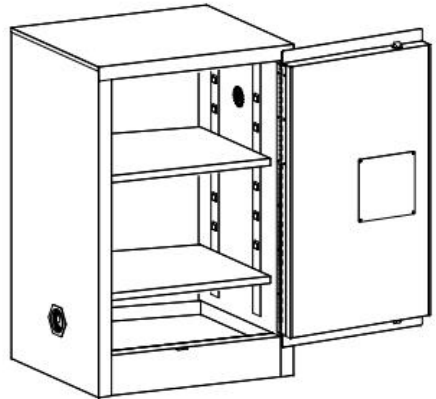
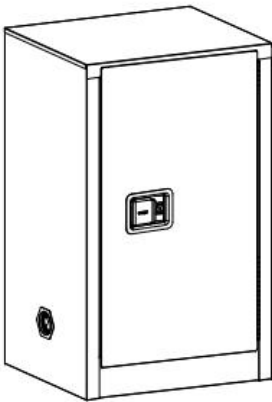
MODEL:SS-13GAL

VEVOR

Affordable. Reliable. Home Improvement.

FLAMMABLE CABINET

MODEL:SS-13GAL

This is the original instruction, please read all manual instructions carefully before operating. VEVOR reserves a clear interpretation of our user manual. The appearance of the product shall be subject to the product you received. Please forgive us that we won't inform you again if there are any technology or software updates on our product.

INSTRUCTIONS

WARNING:

Read this material before using this product. Failure to do so may result in serious injury.

Product Introduction

1. A safety cabinet is a type of storage unit designed to securely store hazardous chemicals, reducing the risk of fire and protecting personnel and equipment.
2. It provides safe storage, containment, and organized management of dangerous chemicals
3. Compact chemical storage cabinets can be placed in your workspace to store flammable substances and organic solvents eliminating the need to go back and forth to a hazardous materials storage room and thus improving efficiency.
4. Even in smaller work spaces, small safety cabinets offer secure storage for flammable liquids and chemicals, reducing raw material waste and saving company costs.

USAGE SAFETY

1. **ATTENTION!** Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in serious injury or even death. The terms "device" or "product" are used in the warning and instructions to refer to Flammable Cabinet.
2. If you are unsure about whether the product is operating correctly or if you find damage, please contact the manufacturer's service centre. Only the manufacturer's service center may make repairs to the product. Do not attempt to make repairs yourself! Keep the device away from sources of fire and heat.

PRODUCT TYPES AND USES

Yellow chemical cabinets are internationally recognized for professional storage of flammable liquid chemicals. These include alcohol, adhesives, gasoline, kerosene, crude oil, ethanol, benzene, toluene, acetone, and more.

PRODUCT FEATURES

1. Complies with OSHA 29 CFR 1910.106 (Occupational Safety and Health Administration regulations) and NFPA CODE 30 (National Fire Protection Association regulations)
2. Constructed with double-layer fire-resistant steel plates, with a 38mm air insulation layer between the layers.
3. High-quality steel plates are point-welded for durability and excellent fire resistance.
4. Three-point interlocking door lock design enhances door sealing
5. Cabinet doors are equipped with dual keys for improved security and regulation of chemical substance management.
6. Professionally standardized warning labels are corrosion-resistant and highly visible.
7. Both sides of the cabinet are equipped with ventilation openings and fire protection devices.
8. Both the interior and exterior of the cabinet are coated with epoxy resin paint.
9. In strict accordance with OSHA regulations, the cabinet body includes electrostatic grounding ports for connecting grounding wires, minimizing the risk of hazards.

INSTRUCTIONS FOR USE

1. Place the explosion-proof safety cabinet on a level surface. If multiple explosion-proof safety cabinets are placed together, the distance between each cabinet should not be less than 15 centimeters.
2. The location for placing the explosion-proof safety cabinet should be away from ignition sources, other heat-generating instruments, and should also be far from splashes of chemicals or slag.
3. At the lower corner of the left exterior wall of the cabinet, there is a static grounding terminal. After properly placing the cabinet, connect the electrostatic grounding wire to discharge any accumulated static electricity from the cabinet.
4. Do not exceed the weight limit of the shelves when placing items on the cabinet shelves. If there are multiple shelves, place larger or heavier items on the bottom shelf and smaller or lighter items on the upper shelf.
5. If the explosion-proof safety cabinet is placed in a confined small space, or if there are strict restrictions on the evaporation of flammable liquids in the cabinet's location, a ventilation system should be connected to the cabinet's vent to eliminate flammable liquids that may accumulate due to evaporation.
6. When placing items inside the cabinet or performing other operations at night or under insufficient light conditions, turn on the surrounding lights or use explosion-proof flashlights for operations.
7. Prohibited actions near the cabinet include lighting candles, alcohol lamps, or using lighters that could potentially cause combustion or explosion.


SAVE THIS MANUAL

PARTS LIST


English Manual*1	
Adjustable foot*4	
Door Key*2	
Ground lead*1	

ASSEMBLY STEP

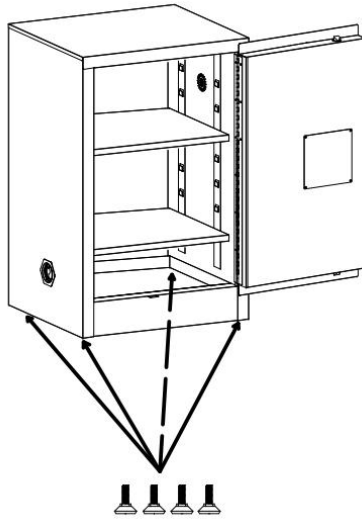When the placement surface is not normal, install the adjusting foot into the reserved thread hole at the bottom, and adjust the height to balance the box.

Adjusting foot installation details:

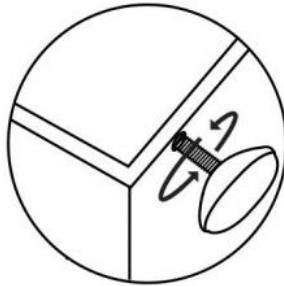PRODUCT PARAMETER

Model	SS-13GAL
Main Color	Yellow
Product Size(W*D*H)	590*460*900mm
Net Weight	44.9kg

* Products such as specifications, appearance, and design are subject to modification without prior notice.

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YH CONSULTING LIMITED.
C/O YH Consulting Limited Office 147, Centurion House, London Road, Staines-upon-Thames, Surrey, TW18 4AX

VEVOR

Affordable. Reliable. Home Improvement.

ARMOIRE INFLAMMABLE

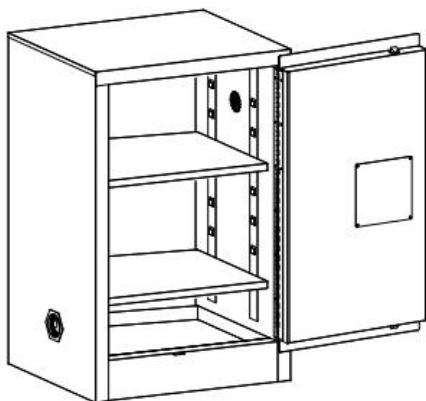
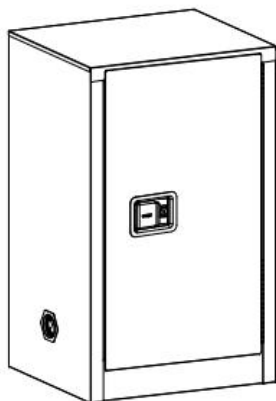
MODÈLE : SS-13GAL

VEVOR

Affordable. Reliable. Home Improvement.

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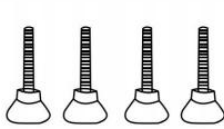
CONSERVEZ CE MANUEL

PARTS LIST

English Manual*1	
Ajustable foot*4	
Door Key*2	
Ground lead*1	

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Importerad till Australien: SIHAO PTY LTD. 1 ROKEVA
Importerad till USA: Sanven Technology Ltd. Suite 250, 9166 Anaheim
E-CrossStu GmbH