# **VEVOR<sup>®</sup>**

**TOUGH TOOLS, HALF PRICE**

Technical Support and E-Warranty Certificate [www.vevor.com/support](http://www.vevor.com/support)

## **LAB DISTILLATION APPARATUS**

### **USER MANUAL**

**MODEL:YHEO-7**

We continue to be committed to provide you tools with competitive price.

"Save Half", "Half Price" or any other similar expressions used by us only represents an estimate of savings you might benefit from buying certain tools with us compared to the major top brands and does not necessarily mean to cover all categories of tools offered by us. You are kindly reminded to verify carefully when you are placing an order with us if you are actually saving half in comparison with the top major brands.

## **SAFETY INSTRUCTIONS**

### **WARNING:**

**Read this material before using this product. Failure to do so can result in serious injury.**

### **Assembly precautions**

1. Assemble only according to these instructions. Improper assembly can create hazards.
2. Wear safety goggles and duty work gloves during assembly.
3. Keep the assembly area clean and well-lit.
4. Keep bystanders out of the area during assembly.
5. Do not assemble when tired or when under the influence of alcohol, drugs or medication.
6. The product capabilities apply to properly and completely assembled products only.
7. Assemble on a flat, level, hard and smooth surface capable of safely supporting the Lab Distillation Apparatus.
8. For additional information regarding the parts listed in the following pages, please refer to the Assembly Diagram of this manual. Unwrap and separate all parts in a clean work area.

### **Use precautions**

1. **DO NOT SIT OR STAND ON THIS ITEM.**
2. This product is not a toy. Do not allow children to play with or near this item.
3. Do not exceed weight capacities.
4. Use only on a flat, level, hard and smooth surface capable of safely supporting a fully loaded Lab Distillation Apparatus.
5. Use as intended only.
6. Inspect before every use; do not use if parts are loose or damaged.

### **SAVE THIS MANUAL**

## INSTRUCTIONS

Thank you very much for choosing this model. Glass fittings should always be handled with care. Please read all of the instructions before using it. The information will help you achieve the best possible results.

Parts name	Number of parts	Description
flask	1	1000ml,24/40 ground glass joint
Two-necked flask	1	250ml,24/40 ground glass joint
Two-necked flask	1	1000ml,24/40 ground glass joint
Three-necked flask	1	500ml,24/40 ground glass joint
Straight condenser	1	/
Coil condenser	1	/
Ball condenser	1	/
Vigreux column	1	/
Vacuum adapter	1	/
Distilling head	1	/
Y-tube	1	/
Thermometer adapter	1	/
Separatory funnel	1	250ml
Glass stopper	5	/
Thermometer	1	200°C
Plastic clip	10	/
Latex tubing	1	2m
PH test paper	2	/

## DIVERSE APPLICATION

Distillation apparatus kit is suitable for chemistry, medical treatment industries, also great for teaching activity, DIY essential oil.

## PRODUCT PARAMETER

<b>Model</b>	<b>YHEO-7</b>
<b>Number of Kits</b>	<b>32</b>
<b>Glass Material</b>	<b>Borosilicate Glass</b>

Manufacturer: Shanghaimuxinmuyeyouxiangongsi  
Address: Shuangchenglu 803nong11hao1602A-1609shi,  
baoshanqu, shanghai 200000 CN.  
Imported to AUS: SIHAO PTY LTD. 1 ROKEVA STREET-  
EASTWOOD NSW 2122 Australia  
Imported to USA: Sanven Technology Ltd. Suite 250,  
9166 Anaheim Place, Rancho Cucamonga, CA 91730

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YH CONSULTING LIMITED.  
C/O YH Consulting Limited Office 147,  
Centurion House, London Road,  
Staines-upon-Thames, Surrey, TW18 4AX

**VEVOR<sup>®</sup>**

**TOUGH TOOLS, HALF PRICE**

**Technical Support and E-Warranty Certificate**

**[www.vevor.com/support](http://www.vevor.com/support)**

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Importerad till AUS: SIHAO PTY LTD. 1 ROKEVA STREET-EASTWOOD
Importerad till USA: Sanven Technology Ltd. Suite 250, 9166
E-CrossStu GmbH