VEVOR[®]

TOUGH TOOLS, HALF PRICE

Technical Support and E-Warranty Certificate www.vevor.com/support

FLAMMABLE CABINET USER MANUAL

We continue to be committed to provide you tools with competitive price.


"Save Half", "Half Price" or any other similar expressions used by us only represents an estimate of savings you might benefit from buying certain tools with us compared to the major top brands and does not necessarily mean to cover all categories of tools offered by us. You are kindly reminded to verify carefully when you are placing an order with us if you are actually Saving Half in comparison with the top major brands.

VEVOR[®]

TOUGH TOOLS, HALF PRICE

FLAMMABLE CABINET

MODEL:SS-889

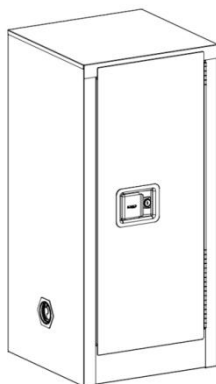NEED HELP? CONTACT US!

Have product questions? Need technical support? Please feel free to contact us:

Technical Support and E-Warranty Certificate
www.vevor.com/support

This is the original instruction, please read all manual instructions carefully before operating. VEVOR reserves a clear interpretation of our user manual. The appearance of the product shall be subject to the product you received. Please forgive us that we won't inform you again if there are any technology or software updates on our product.

Warning-To reduce the risk of injury, user must read instructions manual carefully.

INSTRUCTIONS

Thank you very much for choosing this product. Please read all of the instructions before using it. The information will help you achieve the best possible results.

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual product.

1. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in serious injury or even death. The terms „device“ or „product“ are used in the warning and instructions to refer to Flammable Cabinet.

If you are unsure about whether the product is operating correctly or if you find damage, please contact the manufacturer's service centre. Only the manufacturer's service center may make repairs to the product. Do not attempt to make repairs yourself! Keep the device away from sources of fire and heat.

2. USE GUIDELINES

The product is intended for storing keys out of reach of unwanted persons - it has a key lock. For better organization it has key hangers and an organizer.

CAUTION: Do not store a spare lock key inside the box!

INSTRUCTIONS

1. A safety cabinet is a type of storage unit designed to securely store hazardous chemicals, reducing the risk of fire and protecting personnel and equipment.

2. It provides safe storage, containment, and organized management of dangerous chemicals

3. Compact chemical storage cabinets can be placed in your workspace to store flammable substances and organic solvents eliminating the need to go back and forth to a hazardous materials storage room and thus improving efficiency.

4. Even in smaller work spaces, small safety cabinets offer secure storage for flammable liquids and chemicals, reducing raw material waste and saving company costs.

PRODUCT TYPES AND USES

Yellow chemical cabinets are internationally recognized for professional storage of flammable liquid chemicals. These include alcohol, adhesives, gasoline, kerosene, crude oil, ethanol, benzene, toluene, acetone, and more.

PRODUCT FEATURES

1. Complies with OSHA 29 CFR 1910.106 (Occupational Safety and Health Administration regulations) and NFPA CODE 30 (National Fire Protection Association regulations)
2. Constructed with double-layer fire-resistant steel plates, with a 38mm air insulation layer between the layers.
3. High-quality steel plates are point-welded for durability and excellent fire resistance.
4. Three-point interlocking door lock design enhances door sealing
5. Cabinet doors are equipped with dual keys for improved security and regulation of chemical substance management.
6. Cabinet doors can easily open and close at 180 degrees for convenient and safe operation.
7. 5cm high leak-proof tray prevents accidental spillage of chemical liquids.
8. Professionally standardized warning labels are corrosion-resistant and highly visible.
9. Both sides of the cabinet feature 2-inch diameter vent opening with fire-prevention devices.
10. Unique spill-proof adjustable shelves can be positioned every 6cm, and additional shelves can be added for increased efficiency based on actual chemical storage needs
11. Galvanized steel shelves feature polyethylene spill trays, while the cabinet's bottom has a removable polyethylene liner for easy collection and cleaning of spills (exclusive to blue safety cabinets)
12. Both the interior and exterior of the cabinet are coated with epoxy resin paint.
13. In strict accordance with OSHA regulations, the cabinet body includes

electrostatic grounding ports for connecting grounding wires, minimizing the risk of hazards

INSTRUCTIONS FOR USE

1. Place the explosion-proof safety cabinet on a level surface. If multiple explosion-proof safety cabinets are placed together, the distance between each cabinet should not be less than 15 centimeters.
2. The location for placing the explosion-proof safety cabinet should be away from ignition sources, other heat-generating instruments, and should also be far from splashes of chemicals or slag.
3. At the lower corner of the left exterior wall of the cabinet, there is a static grounding terminal. After properly placing the cabinet, connect the electrostatic grounding wire to discharge any accumulated static electricity from the cabinet.
4. Do not exceed the weight limit of the shelves when placing items on the cabinet shelves. If there are multiple shelves, place larger or heavier items on the bottom shelf and smaller or lighter items on the upper shelf.
5. If the explosion-proof safety cabinet is placed in a confined small space, or if there are strict restrictions on the evaporation of flammable liquids in the cabinet's location, a ventilation system should be connected to the cabinet's vent to eliminate flammable liquids that may accumulate due to evaporation.
6. When placing items inside the cabinet or performing other operations at night or under insufficient light conditions, turn on the surrounding lights or use explosion-proof flashlights for operations.
7. Prohibited actions near the cabinet include lighting candles, alcohol lamps, or using lighters that could potentially cause combustion or explosion.

Section Six: Chemical Storage Methods Store under dry inert gas.

Keep the storage container sealed and place it inside a tightly.

SEALED STORAGE CONTAINER

Store in a cool and dry place.

The storage area must be far from water sources.

The storage area must be far from oxidizers.

Ensure good ventilation and exhaust during working hours.

Must be stored in specialized storage cabinets.



AFTER-SALES SERVICE

The product is covered by a one-year warranty. Please keep the product manual and certific.

TECHNICAL PARAMETERS


Model	SS-889
Colour	Yellow
Size	460*460*900 mm
Material	Q235+Galvanized plate,1.0mm
Door	Single-door,manual

PARTS LIST

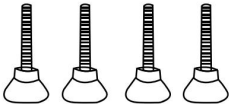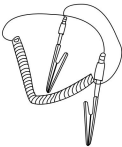
English Manual*1	
Ajustable foot*4	
Door Key*2	
Ground lead*1	

ASSEMBLY STEP

When the placement surface is not normal, install the adjusting foot into the reserved thread hole at the bottom, and adjust the height to balance the box.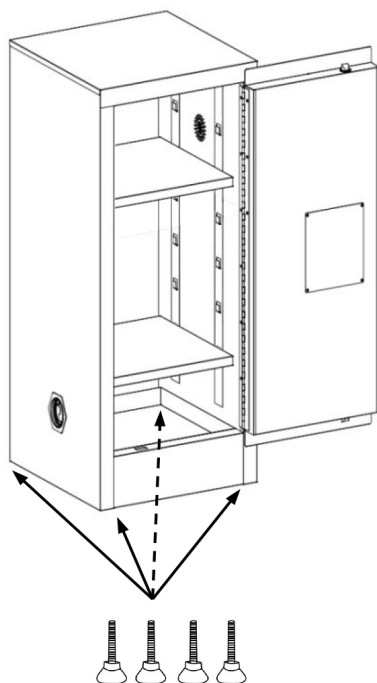

Adjusting foot installation details:

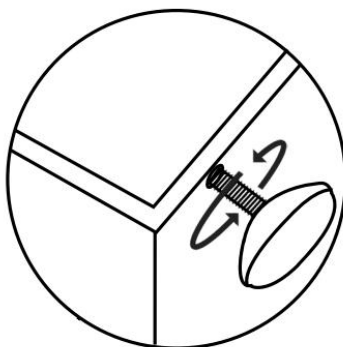Manufacturer: Shanghaimuxinmuyeyouxiangongsi

Address: Shuangchenglu 803nong11hao1602A-1609shi, baoshanqu, shanghai 200000 CN.

Imported to AUS: SIHAO PTY LTD. 1 ROKEVA STREETEASTWOOD NSW 2122 Australia

Imported to USA: Sanven Technology Ltd. Suite 250, 9166 Anaheim Place, Rancho Cucamonga, CA 91730

EC	REP
-----------	------------

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Mainzer Landstr.69, 60329 Frankfurt am Main.

UK	REP
-----------	------------

YH CONSULTING LIMITED.

C/O YH Consulting Limited Office 147, Centurion House, London Road, Staines-upon-Thames, Surrey, TW18 4AX

Technical Support and E-Warranty Certificate
www.vevor.com/support

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Importerad till AUS: SIHAO PTY LTD. 1 ROKEVA STREETEASTWOOD NSW 2122 Australien
E-CrossStu GmbH